



T +44 (0) 2920 854 390
F +44 (0) 2920 854 391

BRUNEL WAY, BAGLAN ENERGY PARK,
NEATH, SA11 2GA, UK
www.ecolabcc.com
infocc@ecolab.com

Klercide™

Filtered Alcohol

Cleanroom filtered 70/30 v/v alcohol, either IPA or Denatured Ethanol, blended with deionised water.

Advantages

■ Easy to use

Convenient range of comfortable and easy to use trigger sprays are available, with lightweight options from 1 to 4 litre. The large droplet size reduces the risk of inhalation.

■ Cost effective

All the sterile alcohol can be dispensed from the bottle so there is no wastage. The SteriShield Delivery System has been validated to give an in-use shelf life of three months, so there is no need to discard unused product.

■ Optimum surface coverage

The fully adjustable trigger spray allows the liquid to be dispensed as either a jet or spray enabling thorough wetting of the surface so that disinfection procedures are effective.



Description

Klercide Filtered Alcohol is available as either 70% v/v Isopropyl Alcohol blended with deionised water or 70% v/v Denatured Ethanol blended with deionised water.

The alcohol blend is 0.2 micron filtered, filled and double bagged in a Grade C (ISO Class 7) cleanroom.

All spray bottles incorporate the patented SteriShield Delivery System and an adjustable trigger spray. The unique way in which the SteriShield Delivery System works, as a closed system, ensures that the integrity of the contents is preserved throughout use.

Available in 1 litre and 4 litre trigger sprays and 5 litre capped bottles.

Protocol

Ready to use so can be sprayed directly onto the surface to be sanitised. Always ensure the whole surface is in contact with the liquid. For optimum results wipe the surface to physically remove any contaminants.

Use biocides safely. Always read the label and product information before use.

Validation

Manufactured by a ISO 9001:2008, ISO 13485:2012 and European Medical Device Directive 93/42/EEC accredited company.

EN test data available on request.

Certificates of analysis with every batch.

2 year shelf life.

Full supporting documentation is available on this product.

Product Codes

Klercide 70/30 IPA

CODE	OLD CODE	LANGUAGES ON LABEL	DESCRIPTION	WATER QUALITY	STERILE / FILTERED	SIZE	UNIT OF SALE
3078430	3036360	GB;DE;FR;IT;ES	Klercide 70/30 IPA	DI	Filtered	1.0 litre	6 bottles x 1.0L
3078410	3055640	GB;SE;NO;FI;PL;DK	Klercide 70/30 IPA	DI	Filtered	1.0 litre	6 bottles x 1.0L
3078400	3036370	GB;DE;FR;IT;ES;NL	Klercide 70/30 IPA	DI	Filtered	4.0 litre	4 bottles x 4.0L
3078420	3036380	GB;DE;FR;IT;ES;NL	Klercide 70/30 IPA	DI	Filtered	5.0 litre	4 bottles x 5.0L

Klercide 70/30 Denatured Ethanol

CODE	OLD CODE	LANGUAGES ON LABEL	DESCRIPTION	WATER QUALITY	STERILE / FILTERED	SIZE	UNIT OF SALE
3078200	3036390	GB;DE;FR;IT;ES;NL	Klercide 70/30 Denatured Ethanol	DI	Filtered	1.0 litre	6 bottles x 1.0L
3078210	3055650	GB;SE;NO;FI;PL;DK	Klercide 70/30 Denatured Ethanol	DI	Filtered	1.0 litre	6 bottles x 1.0L
3038230	3038230	GB;DE;FR;IT;ES;NL	Klercide 70/30 Denatured Ethanol	DI	Filtered	4.0 litre	4 bottles x 4.0L
3078190	3038240	GB;DE;FR;IT;ES;NL	Klercide 70/30 Denatured Ethanol	DI	Filtered	5.0 litre	4 bottles x 5.0L
3078220	3055660	GB;SE;NO;FI;PL;DK	Klercide 70/30 Denatured Ethanol	DI	Filtered	5.0 litre	4 bottles x 5.0L

For more information please contact info@ecolab.com or visit www.ecolabcc.com